

U-SMPS SYSTEM: Sizing and Counting of Nano Scaled Particles

Particles smaller than 0.1 µm have a significant impact not only on health, but also on a wide range of environmental processes such as particulate matter and cloud formation.

Their formation and their effects have long been the focus of scientific attention. For this purpose, systems of different technologies are used for direct or indirect measurement of particle size and concentration.

The **U-SMPS** System manufactured by Palas in Germany meets the diverse normative and technological requirements and have been used successfully for many years in aerosol research and by official measurement networks or private companies.



Application Examples













Principle of Operation

The classification of particles by means of the so-called "electric mobility" and subsequent concentration measurement is a common method for determining the size resolution and distribution of nanoscaled particles in many application areas such as the analysis of synthetically produced nanoparticles or measurement of ultrafine particles of the ambient air.

The determination of size distribution and number concentration is performed by a multi-step procedure:

- Generation of an electrically neutral aerosol (neutralizer):
 For example, XRC 049 (X-ray source), KR-85-370 (radioactive source)
- Continuous, logarithmically ascending size selection (classifier): DEMC SYSTEM, classifier column
- Parallel measurement of raw concentrations (counter):
 UF-CPC SYSTEM, ENVI-CPC SYSTEM, CHARME® SYSTEM
- Raw data inversion (back calculation to original concentrations)
- Diffusion correction (compensation of particle losses due to diffusion processes)



Datasheet

U-SMPS System

All systems provided by Palas for this purpose can be combined and interchanged with each other, thus enabling exact adaptation to the respective application and customer requirements. The open architecture and the interfaces allow third-party systems to be connected in parallel or exclusively to measure the number concentration for comparison.

For example







DEMC 2000 XRC 049 UF-CPC 50

U-SMPS 2050 Measurement range: 4-1,200 nm, $C_N = 0-10^8 \text{ particles/cm}^3$

Standard Combinations

U- SMPS System			1050	2050	2050	2100	1050 X	1100 X	2050 X	2100 X	1700	2700
Classifier	DEMS System	-	х	Х	Х	Х					Х	Х
		Х					Х	Х	Х	Х		
	Columns	1000	х	х			Х	Х			Х	
		2000			Х	Х			Х	Х		Х
Neutralizer	non-radioactive	XRC 049					*	*	*	*		
	radioactive	KR-85									Х	Х
		Ni-63	х	Х	Х	Х						
Counters	Charme	electro- meter									х	х
	ENVI-CPC Sytem	50	х		Х		Х		Х			
		100		Х		Х		Х		Х		
	UF-CPC-System	50	X ¹		X ¹		X ¹		X ¹			
		100		X ¹		X ¹		X ¹		X ¹		

^{*} X Version with integrated XRC049

ENVI-CPC System for continuous ambient air monitoring with UFP-Probe Inlet System integration (EN 16976 / CEN/TS 17434)

UF-CPC System for scientific laboratory measurement (discontinuous/short-duration)

U-SMPS X700 - only for high concentrations

x¹ UF-CPC System as an alternative to ENVI-CPC System

Special Advantages and Benefits

FLEXIBILITY

- Control unit accepts DMAs and counters from different well known manufacturers
- Supports multiple interfaces and remote access

LATEST TECHNOLOGY

- Continuous and fast-scanning principle of measurement
- User-friendly 7" touchscreen with GUI
- Direct graphical visualization of measured data
- Integrated data logger with high storage volume

BEST PRICE-PERFORMANCE RATIO

- Low maintenance
- Reliable functions
- Low operating costs



Palas is a leading developer and manufacturer of high precision instruments for the generation, measurement and characterization of particles in air.

With more than 30 active patents, Palas develops technologically leading and certified fine dust and nanoparticle analyzers, aerosol spectrometers, generators and sensors as well as related systems and software solutions. Palas was founded in 1983 and employs more than 100 people.

Palas GmbH

Siemensallee 84 | Building 7330 | 76187 Karlsruhe Phone: +49 721 96213-0

www.palas.de